

# PTI Technical Note

## In-Situ Calibration on Intracellular $[Ca^{++}]_i$ , $[Na^+]_i$ and pH<sub>i</sub>

Karoly Csatorday<sup>1</sup>, Ram V. Sharma and Ramesh C. Bhalla

<sup>1</sup>Photon Technology International  
University of Iowa, College of Medicine, Iowa City, IA 52242

An expansion of the poster presented at the **THIRD INTERNATIONAL SYMPOSIUM ON CALCIUM AND GENE HYPERTENSION**  
Portland, OR, July 10-14, 1991

### Calcuim

The fluorescence excitation spectrum of the calcium sensitive dye Fura-2 (Grynkiewicz et al. (1985) J. Biol. Chem. 260, 3340-3450) shifts to the shorter wavelengths upon chelating  $Ca^{++}$  ions.

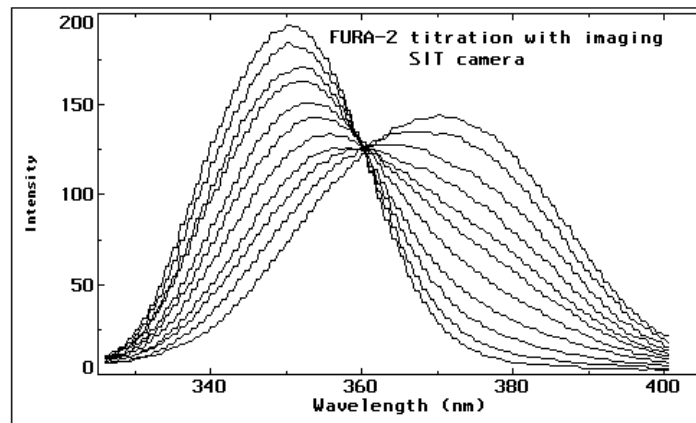


Figure 1: Fluorescence excitation spectra were obtained using a Nikon Diaphot inverted microscope. The UV transmission characteristics determine the peak position of the short wavelength form. In this case the peak is at 348 nm.

The ratio of intensities at 340 and 380 nm reflects the equilibrium between the Ca-bound and Ca-free Fura-2 molecules and can be expressed as  $Ca^{++}$  ion concentration via the analytical expression

$$[Ca^{++}] = K_d \times b \times (R - R_{min}) / (R_{max} - R)$$

where  $k_d$  is the dissociation constant and  $R_{min}$  and  $R_{max}$  are the 340/380 ratios of the Ca-free and Ca-bound forms respectively.  $b$  is the ratio of fluorescence excitation intensities of the two latter forms at 380 nm.  $K_d$  has to be determined independently where-as the other variables are readily accessible from the measurement.

Alternately, by measuring the 340/380 ratios as a function of known free  $Ca^{++}$  concentration, a calibration function called a look-up table (LUT) may be constructed.

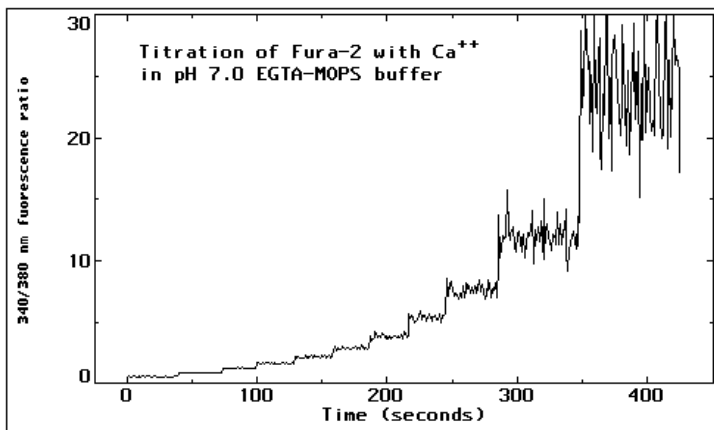


Figure 2: The measured 340/380 fluorescence ratio trace obtained by stepwise titration of Fura-2 with calcium is mapped into concentration space using the analytical formula and the calibration look-up table.

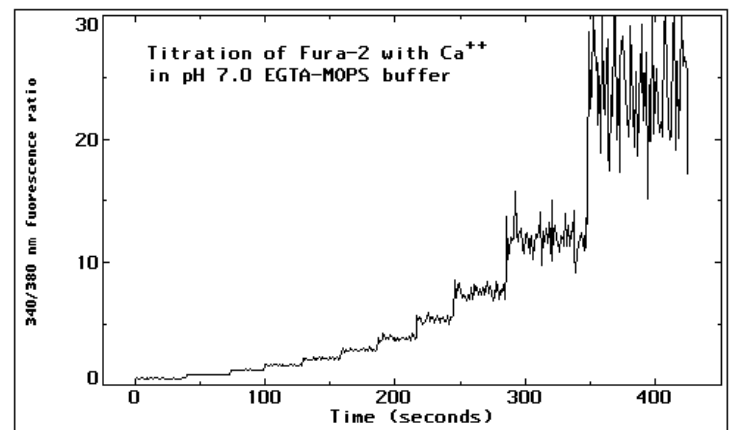


Figure 3: The two traces have been shifted for easy comparison. Congruence of the two traces attests to the viability of the calibration technique via (LUTs).

# PTI Technical Note

## In-Situ Calibration on Intracellular $[Ca^{++}]_i$ , $[Na^+]_i$ and pH<sub>i</sub>

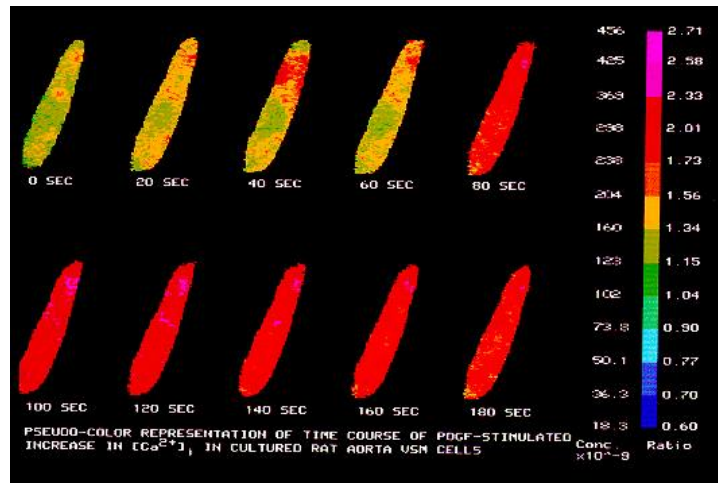


Figure 4: Ratioing fluorescence intensities can be performed on fluorescence images as well. Cells loaded with fluorescent dye are illuminated by alternating 340 and 380 nm light and the resultant fluorescent images are captured by a sensitive video camera and ratioed pixel to pixel. The ratio image is displayed subsequently in pseudo-color with the ratio scale mapped to calcium concentration. (This figure Copyright © 1993 The American Physiological Society. Used by permission.)

### SBFI

A number of researchers attempting to study intracellular  $Na^+$  transients have faced difficulties using the sodium-sensitive dye SBFI. In a study of its behavior in rat aortic vascular smooth muscle cells we reproduced and confirm the results on its spectroscopy in Fluorescence Ratio Imaging of Free  $Na^+$  in Individual Cells by Harootunian A.T. Kao J.P.Y. Eckert B.K. and Tsien R.Y., J. Biol. Chem. 1989, 264, 19458-19467. The imaging hardware of the PTI ImageMaster™ allowed the acquisition of fluorescence excitation spectra directly, using a sensitive video camera that was also used to image the samples.



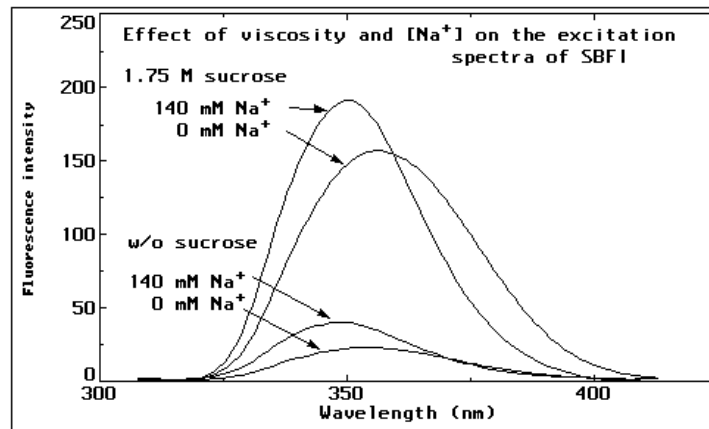


Figure 5: The quantum yield of SBFI fluorescence is 4-6 fold greater in the presence of 1.75 M sucrose than in sucrose-free buffer.

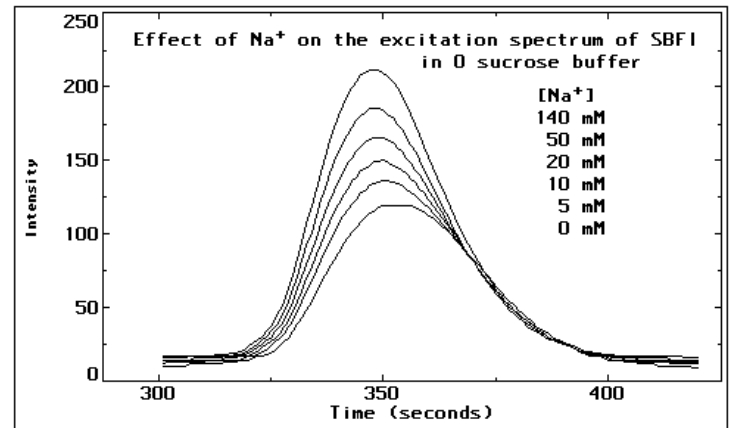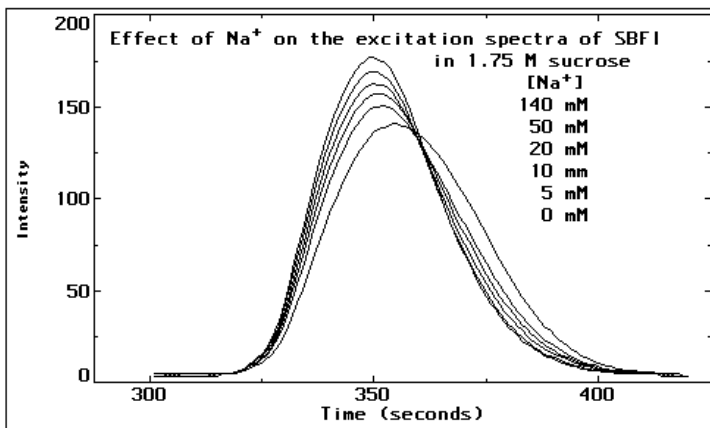

Figures 6 and 7: The spectral shift that occurs upon  $Na^+$  binding is greater in 1.75 M sucrose than in a sucrose-free buffer, facilitating the ratiometric measurement of  $[Na^+]_i$ .

# PTI Technical Note

## In-Situ Calibration on Intracellular $[Ca^{++}]_i$ , $[Na^+]_i$ and pH<sub>i</sub>

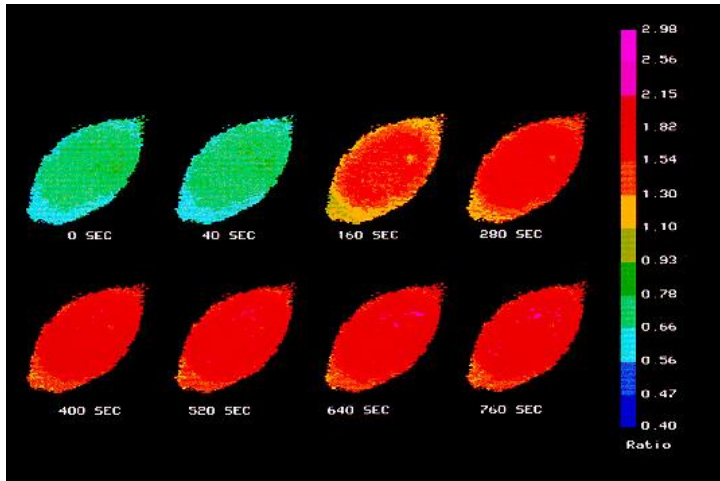


Figure 8: SBFI was observed to indicate higher Na ratios in the nuclear region of VSM cells than in the cytoplasmic region. The effect persisted even in the presence of a mixture of gramicidin/monensin/nigericin which should abolish  $Na^+$ ,  $K^+$  and pH gradients. This leads to the conclusion that the two distinct regions should be calibrated using two different calibration functions, one for the nuclear and another for the cytoplasmic region.

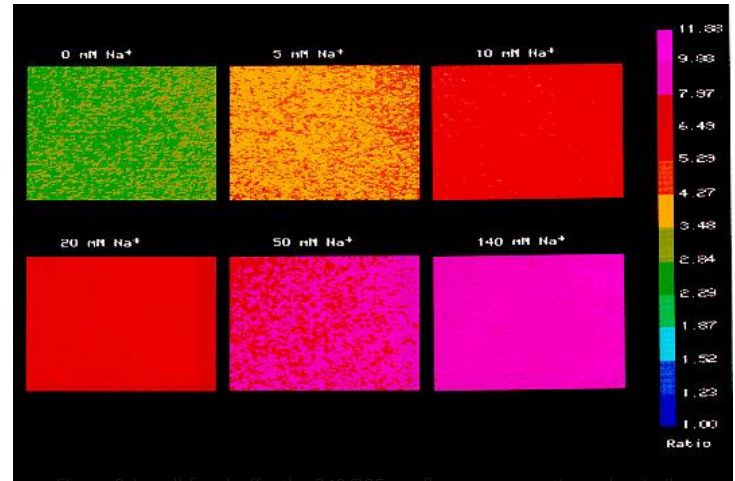


Figure 9: In cell-free buffer the 340/385 nm fluorescence ratio under similar conditions of  $[Na^+]_i$  underwent a much greater change than within VSM cells. Thus such a buffer-based calibration which, if carefully done, can be used in Fura-2 experiments, but will not be adequate to successfully approximate intracellular  $[Na^+]_i$  concentrations.

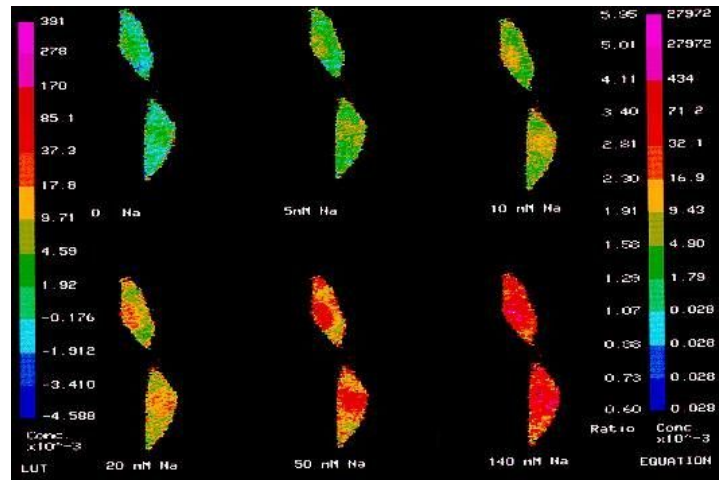


Figure 10: For this composite image, rat aortic vascular smooth muscle cells were clamped at calibrating  $Na^+$  concentrations using gramicidin, monensin and nigericin. The ratio scale was calibrated using a LUT constructed on the basis of Region of Interest photometry from the nuclear region (left scale) as well as based on the concentration equation (right scale).

$R_{max}$  and  $R_{min}$  were 4.3 and 1.1 respectively and  $b$  was measured from the fluorescence excitation spectra to be 1.32. A match in the two scales was achieved in the physiologically meaningful region using a  $K_d$  of 22 mM.

Note: Increasing the number of calibrating points in the look-up table would provide an even closer match in the extreme regions of the scale.

# PTI Technical Note

## In-Situ Calibration on Intracellular $[Ca^{++}]_i$ , $[Na^+]_i$ and pH<sub>i</sub>

### pH

Not having to worry about properly cooling a reflector allows you to concentrate on your work. And, there is no chance of a catastrophic failure ruining an experiment or damaging equipment.

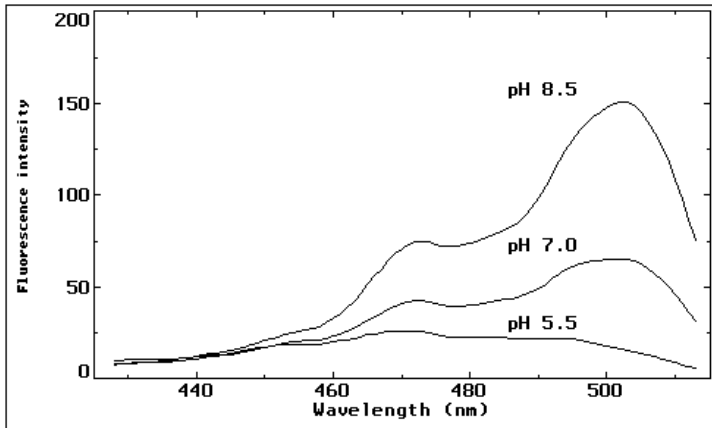


Figure 11: BCECF-loaded VSM cells clamped at calibrating pH values with 10  $\mu$ M nigericin to abolish transmembrane  $[H^+]_i$  gradients were imaged using the technique of fluorescence video microscopy. With the video camera serving as a fluorescence detector, the fluorescence excitation spectra of the pH-sensitive dye BCECF within the cells may be acquired by scanning the excitation monochromators of the ImageMaster™.

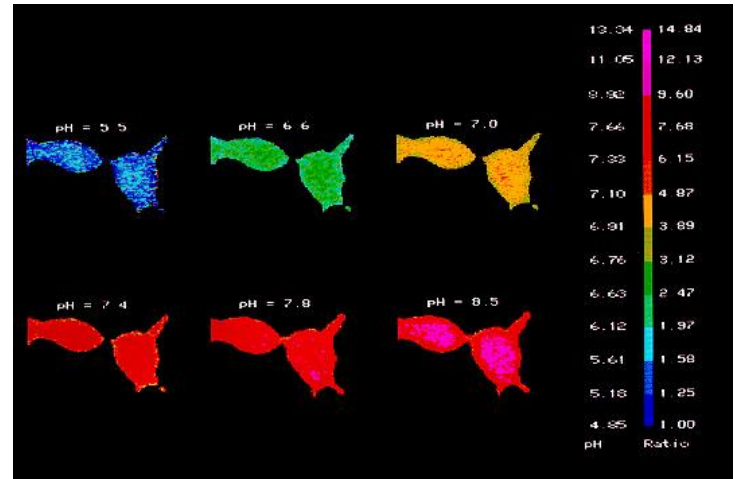


Figure 12: Fluorescence excitation spectra of BCECF at three different pH values were obtained with a SIT video camera. Ratioing the fluorescence images obtained with 500 and 440 nm excitation, a set of pseudocolor images are obtained that serve to calibrate the pH scale.

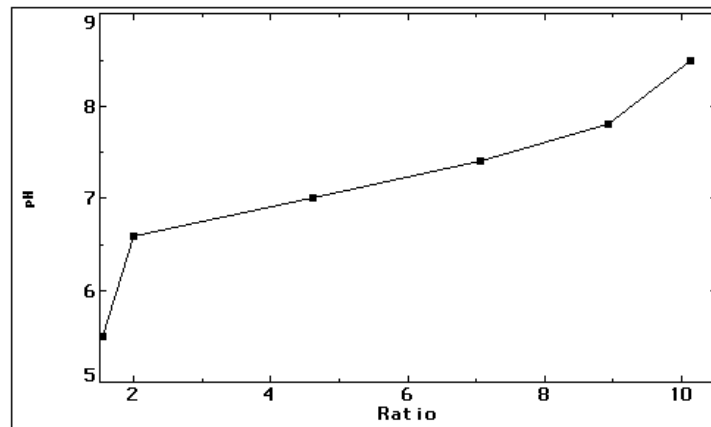


Figure 13: The calibration function is linear in the pH range 6.5-7.5. Changes outside this range may be calibrated by increasing the number of calibrating points in the LUT to span the pH range of interest.

USA: Photon Technology International, Inc., 300 Birmingham Road, PO Box 272 Birmingham, NJ 08011  
Tel: 609-894-4420, Fax: 609-894-1579, E-mail: [marketing@pti-nj.com](mailto:marketing@pti-nj.com), [www.pti-nj.com](http://www.pti-nj.com)

Canada: Photon Technology International, Inc., 347 Consortium Court, London, Ontario, N6E 2S8  
Tel: 519-668-6920, Fax: 519-668-8437, E-mail: [sales@pti-can.com](mailto:sales@pti-can.com)

UK: Photon Technology International, Inc., Unit M1 Rudford Industrial Estate, Ford Road, Ford, West Sussex BN180BF  
Tel: +44 (0) 1903 719555, Fax: +44 (0) 1903 725722. E-mail: [sales@pti-uk.co.uk](mailto:sales@pti-uk.co.uk)

Germany: PhotoMed GmbH, Inninger Str. 1, 82229 Seefeld, Germany,  
Tel: +49 (0) 81 5299 3090, Fax: +49 (0) 81 5299 3098, E-mail: [sales@photomed.com](mailto:sales@photomed.com)



Photon Technology International